

# Bacterial Plating Agar (30 grams)

Catalog #612

EXCITE | EXPLORE | ENGAGE



## Storage: Room Temperature

Most bacterial mediums are prepared at 1.5% agar.

- Media can be prepared by dissolving the nutrient ingredients in water and adjusting the final volume to the desired amount.
- Add the correct amount of agar. Cover the vessel with foil and melt the media on a hot plate or autoclave @121°C for 15-20 minutes.
- Allow media to cool to 55 - 60°C before adding antibiotics or other temperature sensitive additives.
- Swirl media to mix and pour plates. Be careful to avoid contaminants.

1.800.EDVOTEK | [www.edvotek.com](http://www.edvotek.com) | [info@edvotek.com](mailto:info@edvotek.com)