


λ DNA digested with *EcoRI*

Catalog #709

EXCITE | EXPLORE | ENGAGE

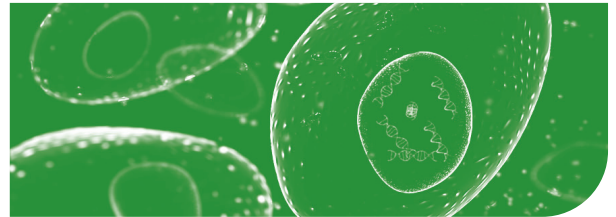

λ DNA digested with *EcoRI*

for 20 gels

The sizes of the λ / *EcoRI* DNA fragments, in base pairs, are 21226, 7421, 5804, 5643, 4878, and 3530. A sample volume of 20 μ l is recommended for ethidium bromide or SYBR Safe staining and 35-40 μ l is optimal for methylene blue staining. Shortly before loading, heat the fragments at 65°C for 3 minutes and quickly chill on ice. The schematic below depicts the λ / *EcoRI* DNA fragments (lane 2) run in parallel with a Standard DNA Gel Marker (cat. #750, lane 1) which is λ / Hind III plus supplemental fragments.

Storage:

Refrigerate or Freeze

